

# User Instructions

81-0201-02 Rev. A



UVP, Inc. 2066 W. 11th Street  
Upland, CA 91786  
(800) 452-6788 / (909) 946-3197  
Fax: (909) 946-3597

Internet: [www.uvp.com](http://www.uvp.com)

Ultra-Violet Product Ltd., Trinity Hall Farm Estate  
Unit 1, Nuffield Rd, Cambridge CB4 1TG UK  
Tel: +44(0)1223-420022  
Fax: +44(0)1223-420561

## Drawer Assembly for the DigiDoc-It System

### Drawer Installation

Place the Benchtop Transilluminator into the drawer. Connect the power cord to the back of the transilluminator and push the plug through the access hole in the back of the drawer. Connect the plug to the power source. Access to the transilluminator switch is via the access port on the front of the drawer.

Place the DigiDoc-it hood enclosure over the drawer. Install the camera. Refer to the DigiDoc-It manual for additional instructions on setting up or using the camera and hood.



*Access to Transilluminator Switch*



*Assemble the System*