

Installation Instructions

81-0210-01 Rev B

Internet: www.uvp.com

UVP, Inc. 2066 W. 11th Street
Upland, CA 91786
(800) 452-6788 / (909) 946-3197
Fax: (909) 946-3597 / E-Mail: uvp@uvp.com



An Analytik Jena Company


Ultra-Violet Product Ltd., Trinity Hall Farm Estate
Unit 1, Nuffield Rd, Cambridge CB1 1TG UK
+44(0)1223-420022 / Fax: +44(0)1223-420561
E-Mail: uvp@uvp.co.uk

Rocker Tray 76-0053-01

Installing the Rocker Tray P/N 76-0053-01

To install the **Rocker Tray** in the HB-1000 or HL-2000, remove the metal bottle carousel from the chamber.

1. Place the tray into the chamber with the coupling arm positioned on the right side.
2. Lift the coupling arm and slide the three pins into the pin hub located on the right side wall of the chamber. Note that the pin lengths are different and can be placed in any position in the pin hub.
3. Slide the tray towards the right so that the pin mount slides against the hub.

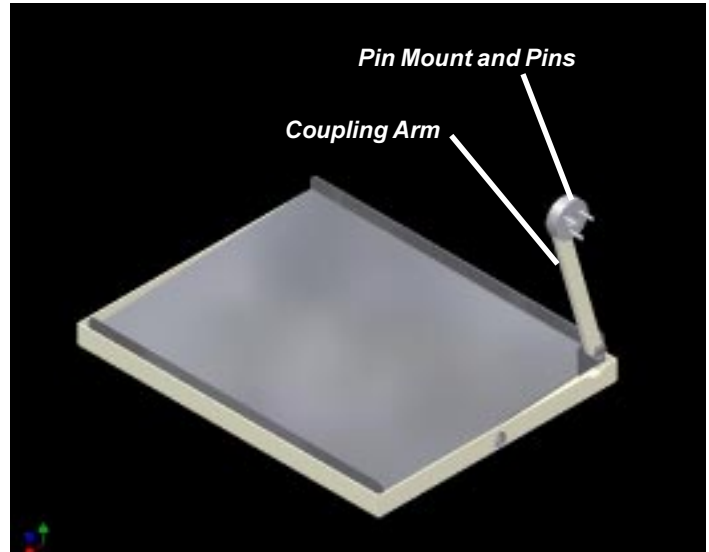

Tray Placement



Hub Location



Pin Mount Placement



Operating the Rocker Tray

After installing the **Rocker Tray**, to operate the tray press the "Motor Control" ON button on the control panel. Adjust the RPM speed from MIN to MAX as required for your experiment.



Control Panel