

# Technical Bulletin

## TB-100 Rev STD

UVP, LLC 2066 W. 11th Street  
Upland, CA 91786  
(800) 452-6788 / (909) 946-3197  
Fax: (909) 946-3597

Web Site: [www.uvp.com](http://www.uvp.com)




Ultra-Violet Product Ltd., Trinity Hall Farm Estate  
Unit 1, Nuffield Rd, Cambridge CB1 1TG UK  
+44(0)1223-420022  
Fax: +44(0)1223-420561

## Power Switch Replacement Instructions

UVP's darkrooms and transilluminators use the same power switch. Follow these steps to replace the power switch in the darkroom or transilluminator:

1. Unplug the darkroom and/or transilluminator from the power source prior to removing the power switch.
2. To remove the power switch from the unit, use a thin flat head screw driver to pry the switch out from housing. Be careful not to scratch the paint of the darkroom or transilluminator in the process.
3. Once the switch is removed, you will see a number of connector wires plugged into the back of it. Normally there are two black and two white wires, with some exceptions. Pull one wire out of the old switch and put it into the same connector location on the new switch. Continue this process for each wire until all are connect to the new switch.
4. Push the new switch back into its position.

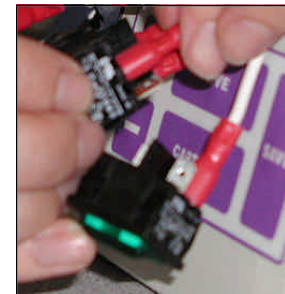

**NOTE: These instructions apply to UVP's Power Switch part numbers:**

53-0126-03	115V
53-0126-01	230V