

UVP White Light Transilluminators (TW-43 and TW-26)

Operating Instructions



Overview

UVP White Light Transilluminators offer the researcher cool white light through an evenly diffused glass surface. The Plexiglas and glass combination is permanently sealed into the unit - double sealed safety fused for durability. Excellent for viewing Coomassie blue stained gels as well as autoradiographs, X-Rays, and microtiter plates. The TW-26 is designed for use with 8 x 10" film sizes; the TW-43 with 14 x 17" film sizes.

Operation

Place the Transilluminator on a level work surface. Be sure that adequate air space exists around the unit. This space allows for proper air circulation through the unit. Plug the female end of the power cord into the Transilluminator. For 230 volt models, or those requiring special power-cord connectors, ensure that the proper configuration of male connector or plug is properly connected to the power cord. Plug the male end of the power cord into a properly grounded electrical outlet. The proper voltage of the Transilluminator is found on the product information label.

Place sample on the filter area. Press the ON/OFF switch to ON. The white lights should be glowing beneath the filter after a momentary flickering during the start period. After viewing/ photographing your sample, turn the unit off.

Cleaning the Transilluminator

The painted surface and filter areas should be cleaned with a damp soft cloth or sponge.

Never use abrasive cleaners as this will damage the filter surface.

ALWAYS DISCONNECT THE TRANSILLUMINATOR FROM ELECTRICAL POWER PRIOR TO CLEANING

Replacing the Tubes

The TW-26 contains two 8-watt tubes; the TW-43 contains four 14-watt tubes. To replace white light tubes:

1. Disconnect the Transilluminator from electrical supply.
2. Remove filter cover by unscrewing retainer screws.
3. Carefully twist the tubes from the lamp holder.
4. Fit the replacement tube into the socket and rotate the tube 1/4 turn until it locks into place.

Replacement Parts

	For TW-26	For TW-43
UVP Filter Assembly	38-0161-04	38-0165-01

UVP Tube, White, 8 Watt 34-0056-01 x2

UVP Tube, White, 14 Watt

34-0059-01 x4

Warranty

Analytik Jena's quality transilluminators are guaranteed to be free of defects in materials, workmanship and manufacture for two (2) years from the date of purchase. Consumable and disposable parts including, but not limited to bottles, tubes and filters, are guaranteed to be free from defects in manufacture and materials for ninety (90) days from date of purchase. If equipment failure or malfunction occurs during the warranty period, AJ shall examine the inoperative equipment and have the option of repairing or replacing any part(s) which, in the judgment of AJ, were originally defective or became so under conditions of normal usage and service.

No warranty shall apply to any instrument, or part thereof, that has been subject to accident, negligence, alteration, abuse or misuse by the end-user. Moreover, AJ makes no warranties whatsoever with respect to parts not supplied by AJ or that have been installed, used and/or serviced other than in strict compliance with the instructions appearing in the operational manual supplied to the end-user.

In no event shall AJ be responsible to the end-user for any incidental or consequential damages, whether foreseeable or not, including, but not limited to property damage, inability to use equipment, lost business, lost profits, or inconvenience arising out of or connected with the use of instruments produced by AJ. Nor is AJ liable or responsible for any personal injuries occurring as a result of the use, installation and/or servicing of equipment. This warranty does not supersede any statutory rights that may be available in certain countries.

Technical Support

AJ offers free lifetime technical support on all of its products and software. Should you have any questions regarding the product's use, operation or repair, contact AJ's offices at the locations below, or visit www.us.analytik-jena.com.

NOTE: A Returned Goods Authorization (RGA) number must be obtained from AJ Customer Service before returning any product.

If you are in North America, South America, East Asia or Australia:	If you are in Europe, Africa, the Middle East or Western Asia:
Call (909) 946-3197 , and ask for Customer Service during regular business days, between 7:00 am and 5:00 pm, PST.	Call +44(0) 1223-420022 , and ask for Customer Service during regular business days between 9:00 am and 5:30 pm.
E-mail your message to: info@us.analytik-jena.com	E-mail your message to: uvp@uvp.co.uk
Fax Customer Service, and send it to (909) 946-3597	Fax Customer Service, and send it to: +44(0) 1223-420561
Write to: Analytik Jena US 2066 W. 11 th Street, Upland, CA 91786 USA	Write to: Ultra-Violet Products Ltd Unit 1, Trinity Hall Farm Estate, Nuffield Road, Cambridge CB4 1TG UK

analytikjena

An Endress+Hauser Company

Analytik Jena US
2066 W. 11th Street
Upland, CA 91786
Phone: (909) 946-3197
Fax: (909) 946-3597

Ultra-Violet Products Ltd.
Unit 1, Trinity Hall Farm Estate
Nuffield Road, Cambridge CB4 1TG UK
Phone: +44(0)1223-420022
Fax: +44(0)1223-420561

Web Site: www.us.analytik-jena.com